


Delineation of Functional Treg Subsets Using Cell Surface Marker Expression

Featured clones:

- Clone V46-1954 is a highly specific anti-mouse CD304 (Nrp-1) antibody that does not cross-react with Nrp-2.
- Clone Y23-1185 is a novel anti-mouse CD39 (NTPDase 1) antibody.

Background

Regulatory T cells (T_{regs}) comprise a heterogeneous group of CD4 T cells that modulate immune responses and maintain immune system homeostasis. Naturally occurring T_{regs} were first characterized within the CD4⁺CD25⁺ T cell subpopulation. These cells express the transcription factor Foxp3, which is essential for T_{regs} stability and immunosuppressive functions. A growing number of cell surface markers support the identification of specific T_{regs} subsets including functional subsets. Neuropilin-1 (Nrp-1 or CD304) is one such marker used to distinguish natural T_{regs} from peripheral and induced T_{regs} in murine models. In both humans and mice, ectonucleoside triphosphate diphosphohydrolase 1 (NTDase 1 or CD39) regulates adenosine-triggered immunosuppressive effects in T_{regs} and has become a novel potential immune checkpoint inhibitor target in cancer immunotherapies.

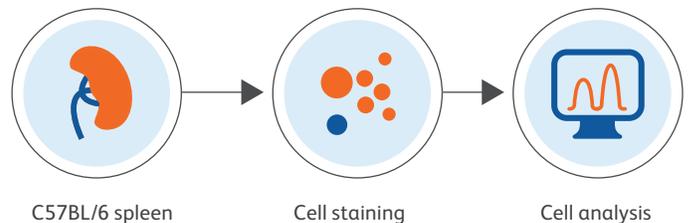

Cell Preparation and Analysis

C57BL/6 mouse spleen was mechanically dissociated and single-cell suspensions were sequentially incubated with Mouse BD Fc Block™ and a cocktail of antibodies. The cells were acquired using a BD FACSymphony™ A5 Cell Analyzer and analyzed in FlowJo™ v10.7.1 Software. Uniform Manifold Approximation and Projection (UMAP) was used for visualization of immune cell populations. Cell populations were annotated based on the expression of lineage-specific markers (Figure 1A). Additionally, CD304 (Nrp-1) expression was shown in a heatmap (Figure 1B). Flow cytometric analysis of the correlated expression of CD304 versus CD25 was derived from CD45⁺CD3⁺CD4⁺ T cells (Figure 2A). Then, CD39 expression was determined within CD304⁻CD25⁻, CD304⁺CD25⁻ and CD304⁺CD25⁺ subsets (Figure 2B).

Laser	Fluorochrome	Specificity	Clone
UV	BUV395	CD8 α	53-6.7
	BUV496	CD45	30-F11
	BUV563	F4/80	T45-2342
	BUV615	CD314	CX5
	BUV661	CD21	7G6
	BUV737	Ly-6C	AL-21
	BUV805	B220	RA3-6B2
Violet	BV421	CD304	V46-1954
	BV510	I-A/I-E	M5/114.15.2
	BV570	CD4	GK1.5
	BV605	TCR γ 6	GL3
	BV650	CD23	B3B4
	BV711	CD27	LG.3A10
	BV750	CD192	475301
Blue	BV786	IgM	R6-60.2
	BB515	CD62L	MEL-14
	BB630	CD11c	HL-3
	BB660	CD357	DTA-1
	PerCP-Cy5.5	IgD	11-26c.2a
	BB755	Ly-6G	1A8
	BB790	NK1.1	PK136
Yellow-Green	PE	CD39	Y23-1185
	PE-CF594	CD93	AA4.1
	PE-Cy5	NKG2A/E/C	20D5
	PE-Cy5.5	CD25	PC61.5
Red	PE-Cy7	CD43	S7
	APC	CD3	145-2C11
	APC-R700	CD19	1D3
	APC-Cy7	CD11b	M1/70

Table 1. Deeper T reg insights from a single tube with broad 29-color immunophenotyping of mouse splenocytes.

BB: BD Horizon Brilliant™ Blue; BUV: BD Horizon Brilliant™ Ultraviolet; BV: BD Horizon Brilliant Violet™

Cat. No	Description	Clone
567112	BD Horizon Brilliant Violet™ 421 Rat Anti-Mouse CD304	V46-1954
567104	BD Pharmingen™ PE Rat Anti-Mouse CD39	Y23-1185

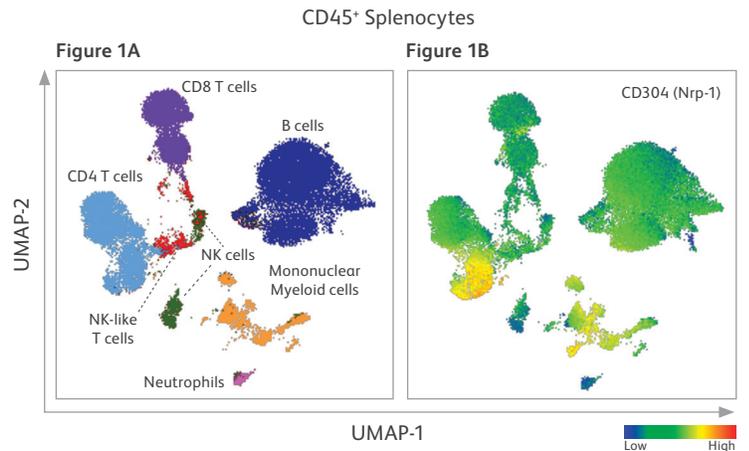


Figure 1. CD304 (Nrp-1) expression most correlated with subset of CD4 T cells and mononuclear myeloid cells.

A. Cell population annotation. B. Heatmap showing CD304 (Nrp-1) expression.

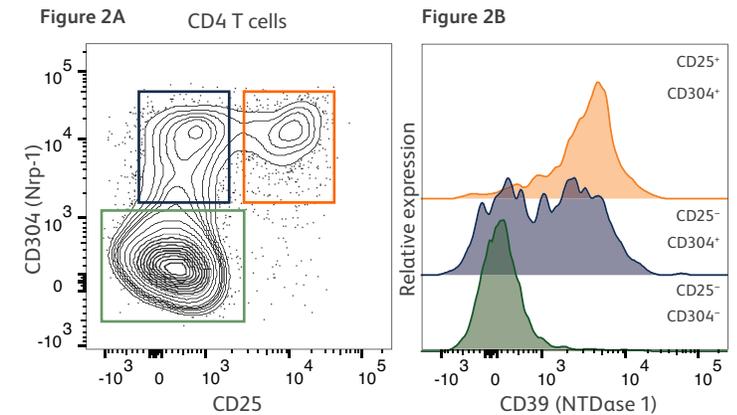


Figure 2. Clean separation of CD4 T cell subsets based on CD304 (Nrp-1) expression and subsequent differential expression of CD39 (NTDase 1).

A. Bivariate contour plots showing CD304 versus CD25 expression on CD45⁺CD3⁺CD4⁺ T cells. B. Histogram overlays showing CD39 (NTDase 1) expression in CD4 T cell subsets.

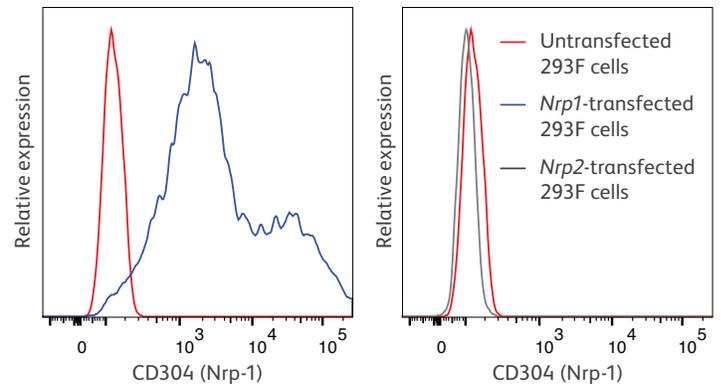


Figure 3. Clone V46-1954 specifically binds CD304 (Nrp-1) and not closely related Nrp-2.

Flow cytometric analysis of CD304 (Nrp-1) expression on transfected cells. 293F cells were left untransfected (red histogram) or transfected with either *Nrp1* (blue histogram) or *Nrp2* (grey histogram). Cells were stained with anti-CD304 (Nrp-1) antibody (clone V46-1954). The histograms show clone V46-1954 is specific to Nrp-1 but not Nrp-2.

BD Life Sciences, San Jose, CA, 95131, USA

bdbiosciences.com

Class 1 Laser Product. For Research Use Only. Not for use in diagnostic or therapeutic procedures.

BD, the BD Logo, FACSymphony, Fc Block, FlowJo, Horizon Brilliant, Horizon Brilliant Violet and Pharmingen are trademarks of Becton, Dickinson and Company or its affiliates. Cy is a trademark of Global Life Sciences Solutions Germany GmbH or an affiliate doing business as Cytiva. CF is a trademark of Biotium, Inc.
© 2021 BD. All rights reserved. BD-30971 (v1.0) 0521

