

Job Aid

BD FACSymphony™ A1 System: Performing daily small particle detector QC

This job aid contains instructions for daily setup of the small particle detector when using the BD FACSymphony™ A1 Flow Cytometer. See the Small particle detector chapter in the *BD FACSymphony™ A1 Flow Cytometer user guide* for more information.

Before you begin

- Create the SPD QC Experiment template described in the BD FACSymphony™ A1 System: Creating the small particle detector QC experiment template job aid.
- Create the SPD setup application settings described in BD FACSymphony™ A1 System: Performing the initial small particle detector setup procedures job aid.
- Start the cytometer and run a performance check.

Preparing the Samples

Preparing the Megamix bead tube:

- Label a clean, capped 12x75 tube *Megamix beads*.
- Mix the Megamix-Plus SSC beads by gently pipetting the mixture.
- Pipet 500 µL of the Megamix-Plus SSC beads into the Megamix beads tube.

Preparing the diluent tube:

- Label a clean, capped 12x75 tube *Diluent*.
- Pipet 3 mL of the 0.2 µm-filtered 1X PBS into the Diluent tube.

Preparing the cleaning tube:


- Label a clean, capped 12x75 tube *Detergent solution*.
- Prepare a tube with 3 mL of a 1.5% dilution of BD® Detergent Solution Concentrate.

NOTE: Never mix BD® Detergent Solution Concentrate and bleach because they create chlorine gas.

Performing small particle detector QC


Setting up the experiment

- ① Verify that your cytometer configuration is appropriate. Your current configuration will need the same as the configuration used to create the SPD QC Experiment template.
- ② Open the SPD QC Experiment template.
 - a. Select a folder in the Browser, then select **Experiment > New Experiment**.
 - b. Select the SPD QC Experiment template and click **OK**.

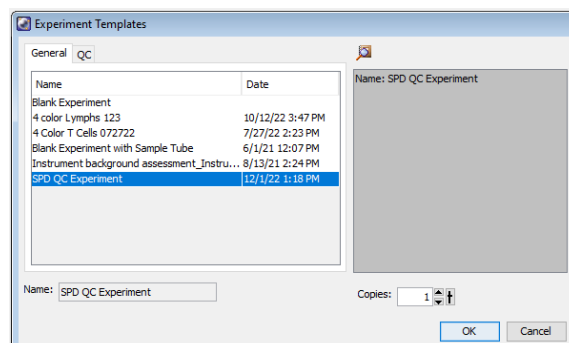

Applying application settings

- ① Apply the SPD setup application settings.
 - a. Right-click the **Cytometer Settings** icon and select **Application Settings > Apply**.
 - b. Select the SPD setup settings you created previously and click **Apply**.
- ② If the Cytometer Settings Mismatch dialog opens, click **Overwrite**.

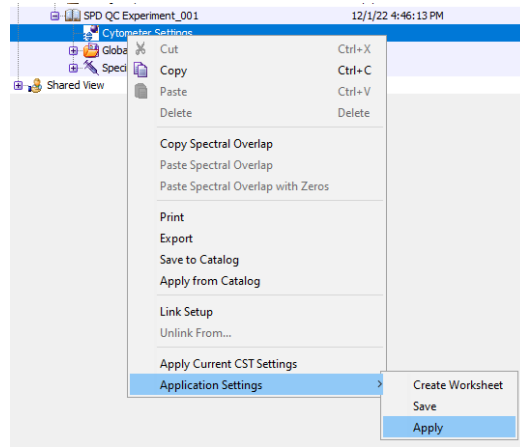


Verifying the application settings

- ① On the cytometer control panel, press the RUN and LOW buttons.

NOTE: For best results when working with small particle detection, the cytometer should stay in run mode with a low flow rate.

- ② Set the current tube pointer to the Megamix beads tube.
- ③ Install the Megamix beads tube onto the cytometer.
 - a. Remove the DI water tube from the SIP.
 - b. Wipe the SIP with a delicate task wipe.
 - c. Install the Megamix beads tube on the SIP. Place the tube support arm under the tube.

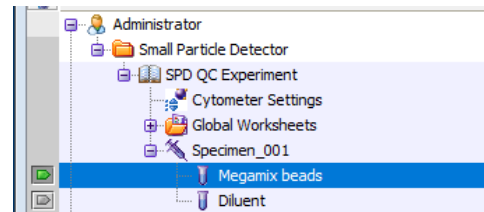

- ④ Click **Acquire Data** in the Acquisition Dashboard.


NOTE: Leave the bead tube on the cytometer until the directions tell you to remove the tube. If you must remove the tube, follow the SPD tube removal procedure.

- ⑤ In the Cytometer window, adjust the SP SSC voltage to place the 200 nm SP SSC-H median at around 100,000, if needed.

- ⑥ Adjust the 200 nm bead gate, if needed.

Verify that the bead peaks are tight and well separated. If not, realign the sample particle detector.

See the BD FACSymphony™ A1 System: Performing the initial small particle detector setup procedures job aid for more information.

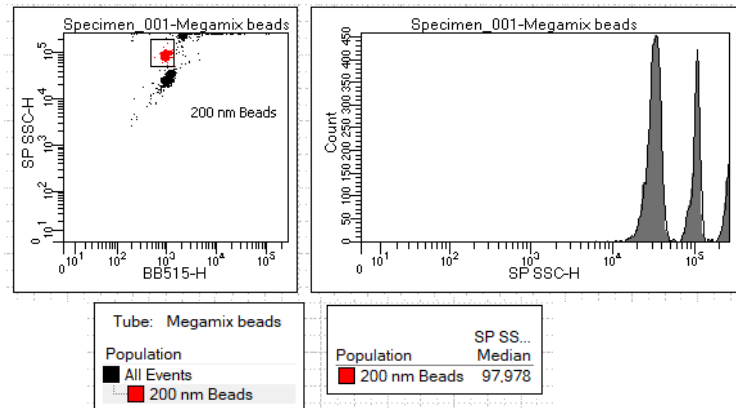


Recording bead data

- ① Click **Record Data**.
- ② When recording is complete, install the Detergent solution tube onto the SIP.
 - a. Remove the Megamix beads tube from the SIP.
 - b. Wipe the SIP with a delicate task wipe.
 - c. Install the Detergent solution tube on the SIP. Place the tube support arm under the tube.
- ③ Run the detergent solution for 15 minutes.

Recording background data

- ① Install the Diluent tube onto the SIP.
 - a. Remove the Detergent Solution tube from the SIP.
 - b. Wipe the SIP with a delicate task wipe.
 - c. Install the Diluent tube on the SIP. Place the tube support arm under the tube.
- ② Set the current tube pointer to the Diluent tube.
- ③ In the Threshold tab, select **SP SSC** as the parameter and verify that the threshold value is set to 200.
- ④ Click **Acquire Data**.
- ⑤ In the Acquisition Dashboard, observe the threshold rate for approximately 5 minutes.
 - The threshold rate should slowly decrease and then plateau.
 - If the threshold rate has not plateaued after 5 minutes, see the Small particle detector troubleshooting section in the BD FACSymphony™ A1 Flow Cytometer user's guide.
- ⑥ Click **Record Data**.
- ⑦ Record for 1 minute, then click **Stop Recording**.

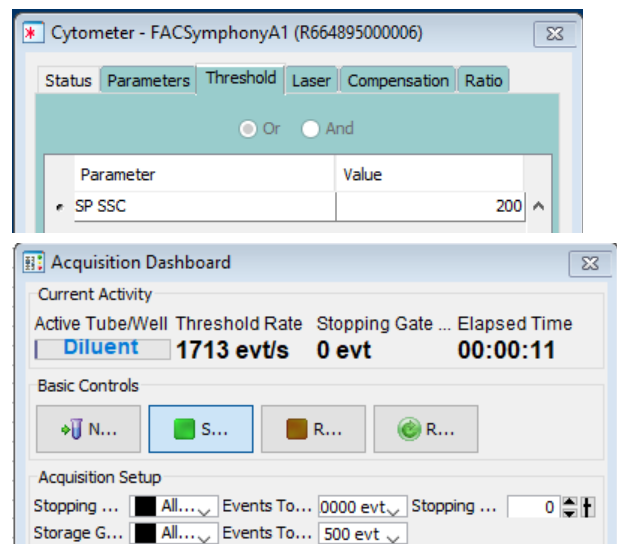

Running your samples

The cytometer is now ready to run your samples.

Tips for running samples:

- Leave the instrument in run mode and low flow rate after starting a small particle experiment, even when waiting between samples.
- Wipe the SIP between samples and avoid leaving the tube support arm to the side for longer than 10 seconds.
- If high background noise is noticed when running samples, clean the fluidics by running a fresh Detergent solution tube for 15 minutes between samples.

This material is for training purposes.
For Research Use Only. Not for use in diagnostic or therapeutic procedures.

BD Life Sciences, San Jose, CA, 95131, USA

bdbiosciences.com

BD, the BD Logo, BD FACSDiva and BD FACSymphony are trademarks of Becton, Dickinson and Company or its affiliates.
©2024 BD. All rights reserved. NPM-1515 (v1.0) 0124

